





Fluoride 100mg/L

DATA SHEET

Instructions for use

- 1. Ensure the pH of the sample roughly 0.5.
- 2. Immerse an indicator strip in the test solution for 1 second.
- 3. Shake off excess solution.
- 4. Wait 60 seconds.
- 5. Compare the colour produced on the strip against the colour chart.

Measuring range

0 - 10 - 25 - 50 - 100mg/L

Application

Fluoride test strips detect fluoride ions in aqueous samples.

Shelf Life

3 years from manufacturing date.

Contents

100 strips

Dimensions

Strip dimension: 83mm x 5mm

Notes

Keep any unused strips in the original container. Do not store where they are liable to be exposed to acid or alkaline gases or vapours. Store in a cool dry place.

E.&O.E.